ON BALANCE™

DZT SERIES

User manual

DZT series USER MANUEL

English

Thank you for purchasing the **DZT series** pocket scale. Please read all operating instructions carefully before use. This electronic scale is a precision instrument. With normal care and proper treatment, it will provide years of reliable service.

- Never load the scale with more than the maximal capacity. Overloading can permanently damage it!
- Avoiding any exposure to extreme heat or cold, your scale works better when operated at normal room temperature. Always allow the scale to acclimate to normal room temperature for at least 1 hour before use.
- Give your scale sufficient warm up time. Usually 30-60 seconds before calibration to give the internal components a chance to stabilize.
- Keep your scale in a clean environment. Dust, dirt, moisture, vibration, air movement and a close distance to other electronic equipment can all cause adverse effects on the reliability and accuracy of your scale.
- Handle with care. Gently apply all items to be weighed onto tray top. Although this scale is designed to be quite durable, try to avoid rough treatment as this may permanently damage the internal sensor.
- Avoid shaking, dropping or otherwise shocking the scale. This is a precision instrument and must be handled with extreme care.
- Only operate the scale on a stable, vibration free surface.

I. OPERATION:
Place scale on a horizontal flat surface
1. Press **ON/OFF** Key.
2. Wait until [0.0] or [0.00] is displayed.
3. Put the object(s) on the weighing platform.
4. Using the **MODE** key, you can switch between the weighing units.

II. TARE WEIGHING:

1. Turn on the scale as described above.
2. Place the "tare item" on the platform.
3. Press **TARE** button and wait until zero is displayed.
4. Add the "net-weight-item".

III. CALIBRATION:

1. Press **ON/OFF** shortly, wait until LCD displays '0.0' or '0.00' and then tap **CAL**.[CAL] is displayed.
2. Tap CAL again, required calibration weight will be displayed.
3. Place the requested weight on the center of platform, wait 3-4 seconds. The LCD display reads [PASS]
4. Remove the weight and press **ON/OFF** to turn off the scale.

Finished
Repeat this process if something went wrong.

Please check the availability of weights at the store you bought your scale from.

IV. SPECIFICATIONS:
Capacity: 600g/21.16oz/19.29oz/385.84wt
9259gn/3000.0ct
Readability: 0.1g/0.01oz/0.01oz/0.1dwt
1gn/0.5ct
Units: g/oz/gn/ct/ozt/dwt
Capacity: 100g/3.527oz/3.215oz/64.30dwt
1543.2gn/500.00ct
Readability: 0.01g/0.001oz/0.001oz/0.01dwt
0.1gn/0.5ct
Units: g/oz/gn/ct/ozt/dwt

Auto off time: 180 seconds
Operation temp: 10-30°C
Display: LCD 5 1/2 digits
Power: 2xAAA Batteries
Dimensions: 128x77x19 mm
Weight: 130g

V. TROUBLESHOOTING:

The primary reasons for inaccuracy or malfunction are low batteries, incorrect calibration, overload or operating on an unstable surface. Please keep these in mind and maintain and operate your scale properly. The scale is a precise instrument and must be handled with the utmost care and caution.

Display	Elimination of errors
Nothing, Battery-symbol, [lo] or [8888]	Change the batteries.
[OUT2] [EEEE]	Recalibrate the scale as shown above.
[LLLL]	You are overloading the scale. Remove the excessive load immediately!
[UNST]	The delicate weighing load cells have been damaged by missing the scale (i.e. it was shocked, dropped or overload.) You can try recalibrating the scale. If the load cell has not been damaged too much it will work again after recalibration.
	Use the scale in a more stable position

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